

Am Jägersberg 5-7
24161 Altenholz
Germany

phone: + 49 - 4 31 - 3 69 60 - 0
fax: + 49 - 4 31 - 3 69 60 21
mail: info@hydrobios.de /
sales@hydrobios.de
web: www.hydrobios.de



Apparatebau GmbH

Operating Instruction

Repair Kit for Net Bucket (Cat. No. 438 025-001 and 438055-001)

The repair kit for net buckets contains a mesh in the ordered mesh size (here 780µm), 4 tie raps and 4 stripes of double sided self-adhesive tape.



Additionally you will need a sharp cutter or a scalpel and some sort of cleaning solvent to degrease the surfaces of the net bucket (e.g. acetone or isopropanol).



As a first step, the old mesh has to be removed from the net bucket, Please cut off the old tie raps and remove the mesh from the self-adhesive tape. Afterwards please also remove the old self-adhesive tape from the bucket and clean the bucket thoroughly with the cleaning solvent.



Apply the self-adhesive tape as shown in the two pictures below:



Take of the protection paper of the self-adhesive tape but without removing it completely.




Now apply the mesh, beginning in the middle of the window. Stretch the mesh a bit towards the ends of the bucket in order to avoid wrinkles.




Finally remove the rest of the protection paper and press the rest of the mesh to the self-adhesive tape. Repeat applying pressure along the self-adhesive tape once again.



Now apply the tie raps, beginning in the outer notches. Fix them firmly.

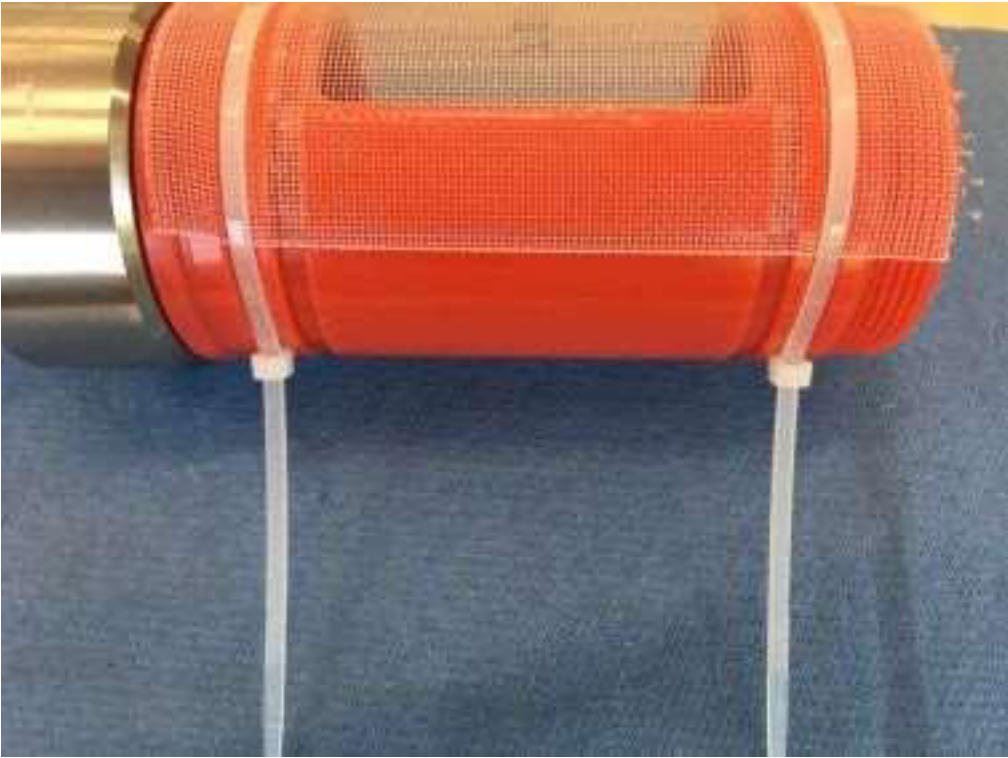

Now install the tie raps to the inner notches and fix them firmly.

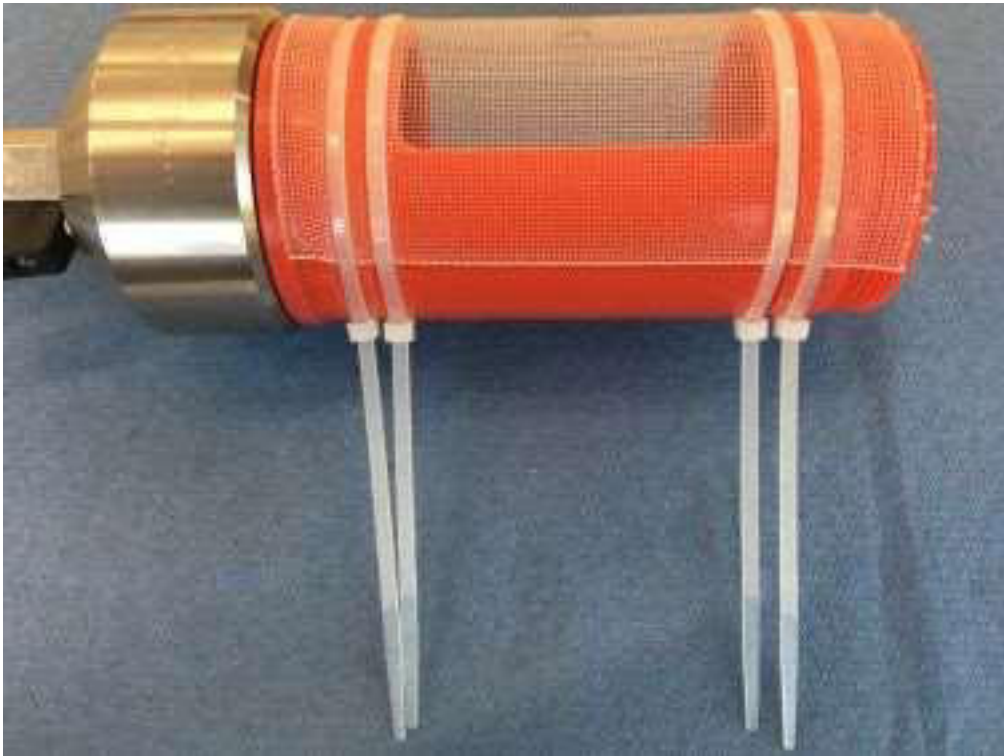

Finally, the overlapping mesh has to be cut-off carefully. In addition, the flaps of the tie raps have to be cut-off, Please mind that the cut should be as close as possible to the lock of the tie raps in order to avoid sharp edges which might damage the net during use.



Revision 1, 30.04.2019